

🐭 H&E staining frozen sections

1. Fixation:
 - Acetone at -20 °C, for 3 min.
 - 80% Methanol at 4°C, for 5 min
2. PBS 1X 10 min.
3. Hematoxylin 1 min (filter before use).
4. Distillated H₂O, for 5 min.
5. Eosine 30 sec.
6. Ethanol 70%, for 5 min.
7. Ethanol 80%, for 5 min.
8. Ethanol 90%, for 5 min.
9. Ethanol 100%, for 5 min.
10. Xylene for 5 min.
11. Xylene for 5 min.
12. Dry completely.
13. Wet slides with Xylene.

NOTE: to mount slides use Permount.

